

56 Sparta Avenue • Newton, New Jersey 07860 (973) 300-3000 Sales • (973) 300-3800 Fax

PA44LE-16 - March 8, 2018

Item # PA44LE-16 was discontinued on March 8, 2018. For informational purposes, this is a copy of the website content at that time and is valid only for the stated product.

# PIEZOELECTRIC RING STACKS AND CHIPS, 2.6 MM TO 9.0 MM TRAVEL



Click to Enlarge

A silver dot indicates the positive electrode (above, left). There is no mark next to the

negative electrode (above, right).

## Hide Overview

### OVERVIEW

# Features

- Ring Cross Section with 8.3 mm
   OD and 3.0 mm ID
- Chips with Bare Electrodes or Pre-Attached Wires
- Discrete Stack with Flat Ceramic End Plates and Pre-Attached Wires
- Drive Voltage Range of 0 150
  V
- For Use in Open-Loop Setups
- Un-Wired Chips in Packs of 16 Available
- Ideal for Vacuum and OEM
- Applications
- · Custom Options Available; Please Contact Tech Support

Thorlabs' offers our piezoelectric ring actuators as a chip with bare electrodes, chip with pre-attached wires, and as a discrete stack consisting of multiple chips bonded via epoxy and glass beads. They have outer diameters of 8.3 mm and inner diameters of 3.0 mm. The wired and unwired chips provide a maximum free stroke of 2.6 µm, and the discrete stack provides 9.0 µm of travel. The central through hole makes these products ideal for laser tuning and micro-dispensing applications. They are available with a drive voltage range of 0 - 150 V; for a complete list of specifications, see the tables below. For convenience, 75 mm wires are attached to the PA44LEW chip and the PK44LA2P2 stack. The flat endfaces of the PK44LA2P2 stack are fitted with ceramic end plates.

#### Circular vs. Square Outer Dimensions

The circular geometry of these ring piezos allow the load to be uniformly distributed across the surface, while piezo actuators with square outer dimensions experience stress concentration at their sharp corners. Cracks, which can lead to failure of the actuator, form more easily in areas of concentrated stress. Greater immunity from stress-induced cracking allows ring chips to be fabricated with thinner

active layers than the square piezoelectric actuators, which results in the ring piezo actuators achieving higher displacements for the same driving voltages. Our design is optimized to provide ring piezo actuators with the longest stroke. For a driving voltage of 150 V, the free stroke per length ratio is approximately 44% greater for ring piezo chips and 28% greater for the ring piezo stacks than for the similar versions with square outer dimensions. Both types possess compact dimensions and light weight, and they can easily be integrated into systems for precision movement. Piezo actuators with square outer dimensions may be preferred over ring piezos when an actuator with a rectangular shape is more easily accommodated, such as in a square corner, or when lower-cost actuators are desired.

#### Piezo Chip Construction and Operation

Piezo chips are fabricated from a series of stacked piezoelectric ceramic layers. An electrode is screen-printed on each layer, and the printed layers are isostatically pressed together to create a chip. The electrodes are electrically in parallel. Through a precision grinding process, the accuracy of the design height is ensured to better than ±5 µm. This high accuracy makes it significantly easier to design devices around our piezoelectric chips and stacks. Each chip has an insulating ceramic layer sealing the outer and inner diameters, as well as the flat top and bottom surfaces. The drive voltage is applied to two screen printed electrodes located on the outer diameter of the ring. For these chips and stacks, a silver dot indicates the positive electrode that should receive positive bias, as is shown in the image at the top of this section. In the case of the wire actuators, the wire soldered to this electrode is red.

Piezo Selection Guide <sup>a</sup>
Piezo Chips
Square
Square with Through Hole
Round
Ring
Shear
Benders
Piezo Stacks
Discrete, Square
Discrete, Square with Through Hole
Discrete, Round
Discrete, Ring
Discrete, Shear (1D to 3D Positioners)
Co-Fired, Square
Co-Fired or Discrete, Square with Strain
Gauges
Piezo Actuators
Mounted

 For more information about the design and function of piezoelectric chips, please see our piezoelectric tutorial. The maximum displacement of these actuators is achieved when they are preloaded with the maximum displacement load, which is specified for each product. The actual value of the maximum displacement varies for each item and must be experimentally determined; however, the maximum displacement will always be larger than the free stroke displacement. Please note that when mounting a load onto the piezoelectric chip, the force should be directed along the actuator's axis of displacement. For more details see the *Operation* tab.

Piezo chips with custom dimensions, voltage ranges, and coatings are available. Additionally, customers can order these piezo chips in high-volume quantities. Please contact Tech Support for more information.

#### Thorlabs' In-House Piezoelectric Manufacturing

Our piezoelectric chips are fabricated in our production facility in China, giving us full control over each step of the manufacturing process. This allows us to economically produce high-quality products, including custom and OEM devices. A glimpse into the fabrication of our piezoelectric chips follows. For more information about our manufacturing process and capabilities, please see our Piezoelectric Capabilities page.

- · Build Blocks from Flexible Sheets of Lead Zirconate Titanate (PZT) Powder
  - Screen Print Electrodes on Each Individual Sheet
  - Layer the Printed Sheets One Top of AnotherConsolidate the Layered Sheets in an Isostatic Press
- · Dice the Block into Individual Elements
- · Purge Solvent and Binder Material Residues by Heat Treating the Elements
- · Sinter the Elements to Fuse the Piezoelectric Pressed Powder and Grow PZT Crystals
- · Lap the Elements to Achieve Tight Dimensional Tolerances: ±5 µm for Each Element
- · Screen Print the Outer Electrodes on the Elements
- · Align the Individual PZT Crystals Along the Same Axis by Poling the Elements

# Hide Operation

### OPERATION

# **Operation Notes**

#### Power Connections

A positive bias should be applied across the device. The positive electrode should receive positive bias, and the other electrode should be connected to ground. Applying a negative bias across the device may cause mechanical failure. For products that ship with wires attached, the positive wire is red, as can be seen in the product images. For pre-wired chips, the red wire is attached to the electrode that is identified by a + mark, as is depicted in Figure 1. (On some devices, a silver dot indicates the positive electrode.) On stacks, the red wire is attached to electrode on the side of the stack indicated by the line of + marks (or silver dots), as is depicted in Figure 2. The wire that should be grounded is black, and it is attached to the electrode on the side of the chip opposite the side with the positive electrode.

#### Preloading

The maximum displacement of these actuators is achieved when they are preloaded with the maximum displacement load, which is specified for each product. The actual value of the maximum displacement varies for each item and must be experimentally determined; however, the maximum displacement will always be larger than the free stroke displacement. Preloading increases the length of the actuator's stroke because the poling process performed during fabrication does not align all ferroelectric grains in the piezoelectric material in the same direction. Preloading the actuator mechanically forces many of the mis-aligned grains into a more ideal alignment. Applying a driving voltage across the piezo material causes the orientations of the ferroelectric grains to rotate so they become aligned with the applied field, and this results in a dimensional change of the piezo material. When more ferroelectric grains are initially aligned in the same direction, the dimensional change of the piezo material in response to the applied driving voltage is greater. Preloads greater than the optimal maximum displacement load result in displacements less than the maximum displacement, as higher loads oppose the switching of the grain orientations in response to the applied driving voltage.

#### Soldering Wire Leads to the Electrodes

If wire leads must be reattached to the electrodes, a soldering temperature no higher than 370 °C (700 °F) should be used, and heat should be applied to each electrode for a maximum of 2 seconds. Solder the lead to the middle of the electrode and keep the region over which heat is applied as small as possible.

#### Interfacing a Piezoelectric Stack with a Load

Piezoceramics are brittle and have low tensile strength. Avoid loading conditions that subject the actuator to lateral, transverse, or bending forces. When applied incorrectly, an external load that may appear to be compressive can, through bending moments, cause high tensile stresses within the piezoelectric device. Improperly mounting a load to the piezoelectric actuator can easily result in internal stresses that will damage the actuator. To avoid this, the piezoelectric actuator should be interfaced with an external load such that the induced force is directed along the actuator's axis of displacement. The load should be centered on and applied uniformly over as much of the actuator's mounting surface as possible. When interfacing the flat surface of a load with an actuator capped with a flat mounting surface, ensure the two surfaces are highly flat and smooth and that there is good parallelism between the two when they are mated. To attach a load to the piezo stack, we recommend using an epoxy that cures at a temperature lower than 80 °C (176 °F), such as our 353NDPK or TS10 epoxies or Loctite<sup>®</sup> Hysol<sup>®</sup> 9340. Loads should be mounted only to the faces of the piezoelectric stack that translate. Mounting a load to a non-translating face may lead to the mechanical failure of the actuator.

#### **Operating Under High-Frequency Dynamic Conditions**

It may be necessary to implement an external temperature-control system to cool the device when it is operated at high frequencies. The maximum operating temperature of these devices is 130 °C (266 °F), and high-frequency operation causes the internal temperature of the piezoelectric device to rise. The dependence of the device temperature on the drive voltage frequency for each product can be accessed by clicking the Info icons,  $\mathbf{0}$ , below. The temperature of the device should not be allowed to exceed its specified maximum operating temperature.



Chips After Binder Burnout and Sintering

÷

Negative Wire (~75 mm Long

Positive Wire

Click to Enlarge Figure 1 : Schematic of TA0505D024W Piezo Chip

Click to Enlarge Figure 2: Schematic of PK4DMP2 Piezo Stack Showing the Line of + Marks Close to the Positive Electrode

owing the + Mark Close to the Positive Electrode



Click to Enlarge Dicing the PZT Block into Individual Elements



#### Estimating the Resonant Frequency for a Given Applied Load

A parameter of significance to many applications is the rate at which the piezoelectric actuator changes its length. This dimensional rate of change depends on a number of factors, including the actuator's resonant frequency, the absolute maximum bandwidth of the driver, the maximum current the piezoelectric device can produce, the capacitance of the piezoelectric stack, and the amplitude of the driving signal. The length of the voltage-induced extension is a function of the amplitude of the applied voltage driving the actuator and the length of the piezoelectric stack. The higher the capacitance, the slower the dimensional change of the actuator.

Quick changes in the applied voltage result in fast dimensional changes to the piezoelectric stack. The magnitude of the applied voltage determines the nominal extension of the stack. Assuming the driving voltage signal resembles a step function, the minimum time,  $T_{min}$ , required for the length of the actuator to transition between its initial and final values is approximately 1/3 the period of resonant frequency. If there is no load applied to the piezoelectric stack, its resonant frequency is  $f_{c}$  and its minimum response time is:

$$T_{min} \cong \frac{1}{3f_o}$$


After reaching this nominal extension, there will follow a damped oscillation in the length of the actuator around this position. Controls can be implemented to mitigate this oscillation, but doing so may slow the response of the actuator.

Applying a load to the actuator will reduce the resonant frequency of the piezoelectric stack. Given the unloaded resonant frequency of the actuator, the mass of the stack, m, and the mass of the load, M, the loaded resonant frequency ( $f_o$ ) may be estimated:

$$f_o' \cong f_o \sqrt{\frac{m_{/3}}{m_{/3} + M}}$$

### Co-Fired Piezo Chips, Co-Fired Piezo Stacks, and Discrete Piezo Stacks

Generally speaking, Thorlabs' co-fired piezo chips and stacks are constructed similarly. In both cases, a structure is built up from alternating electrode layers and green-state lead zirconate titanate (PZT) piezoelectric layers. The assembled structure is then sintered into a single monolithic unit. Supply electrodes of opposite polarity are attached to opposite sides of the structure. Each internal electrodes of opposite board to one or the other supply electrode, such that no two adjacent internal electrodes have the same polarity. The most significant differences between the co-fired chips and stacks arise from the way each internal electrode is electrically isolated from the supply electrodes of opposite polarity. The different electrical insulation approaches influence the mechanical properties of the actuators. Actuators fabricated using the two different approaches are diagrammed in Figure 3.

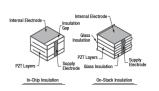

Click to Enlarge Figure 3: Diagram of Piezo Stack Insulation Methods: (a) In-Chip Insulation Used in Standard Chips and Discrete Stacks, (b) On-Stack Insulation Used in Co-Fired Stacks

In the case of the chips (In-Chip Insulation), the internal electrodes of opposite polarities alternate. Each internal electrode layer is shorter than the full width of the piezo layer. All electrodes of one

polarity have edges that are flush with one side of the chip, and all electrodes with the opposite polarity are flush with the opposite side of the chip. Because the electrode does not extend all the way to the opposite edge, the far end of the electrode is completely surrounded by PZT material. The PZT material enclosing the end of the electrode is insulating, which electrically isolates this electrode from the supply electrode of opposite polarity. This approach to electrically insulating the electrode screates a region of stress at the edge of the electrode. The stress arises both due the abrupt change in thickness on either side of the electrode edge, as well as the tensile stress created when the PZT material sandwiched between electrodes responds to an applied voltage drive signal, but the insulating PZT material beyond the edge of the electrodes not. This stress limits the maximum height of chips manufactured using this approach. The height of chips are limited to ensure internal stresses are low and do not affect lifetime or performance. Chips are sealed in a ceramic layer that offers superior resistance to humidity and heat than epoxy resin coatings.

One way of increasing the height, and therefore the maximum stroke, of piezo actuators based on these chips is to fabricate discrete piezo stacks. These are manufactured by bonding multiple chips together in series using a glass-bead epoxy. Discrete stacks can be fabricated to substantially longer lengths than cofired chips or stacks, and this allows them to achieve higher maximum displacements while maintaining sub-millisecond response times and a low drive voltage range. As the constituent chips are sealed within a ceramic barrier layer, discrete stacks have superior resistance to humidity and heat than co-fired stacks, which are sealed in an epoxy resin coating.


In the case of co-fired stacks (On-Stack Insulation), the electrodes extend across the full width of the PZT layers. The edges of the electrodes are flush with all four sides of the stack, including the side with the supply electrode of opposite polarity. The edge of the internal electrode is insulated from that supply electrode by a layer of glass filament applied to the side of the stack. Precision localized application of the glass filament ensures that the electrode edge is electrically isolated from the supply electrode, and that the filament is applied over minimal surface area; the ability of the supply electrode to make electrical connections to the desired internal electrodes is not affected, and the small amount of applied glass filament does not affect the operation of the actuator. With their full-width electrodes, piezo actuators made using this insulating approach are characterized by homogeneous internal stress. Co-fired stacks can therefore be fabricated with greater heights than chips fabricated using the in-chip insulation approach. Co-fired stacks also have a higher percentage of active PZT material than the discrete stacks, which include inactive bonding layers of glass-bead epoxy. They are coated in an epoxy resin.

#### Estimating Device Lifetime for DC Drive Voltage Conditions

The lifetime of a piezoelectric device is a function of the operating temperature, applied voltage, and relative humidity conditions. Lifetimes are reduced as a consequence of humidity-driven electrolytic reactions, which occur at the electrolects of the piezoelectric devices when a DC voltage is applied. These reactions both generate hydrogen and result in metal dendrites growing from the cathode towards the anode. The hydrogen liberated by the electrolytic reaction chemically reacts with and degrades the piezoelectric material. Dendrites that grow to electrically connect the cathode and anode result in increasing levels of leakage current. Failed piezoelectric devices are defined as those that exhibit leakage current levels above an established threshold.

A ceramic moisture-barrier layer that insulates Thorlabs' piezoelectric devices on four sides is effective in minimizing the effects of humidity on device lifetime. As there is interest in estimating the lifetime of piezoelectric devices, Thorlabs conducted environmental testing on our ceramic-insulated, low-voltage, piezoelectric actuators. The resulting data were used to create a simple model that estimates the mean time to failure (MTTF), in hours, when the operating conditions of humidity, temperature, and applied voltage are known. The estimated MTTF is calculated by multiplying together three factors that correspond, respectively, to the operational temperature, relative humidity, and fractional voltage of the device. The fractional voltage is calculated by dividing the operational voltage by the maximum specified drive voltage for the device. The factors for each parameter can be read from the following plots, or they may be calculated by downloading the plotted data values and interpolating as appropriate.

In the following trio of plots, the solid-line segment of each curve represents the range of conditions over which Thorlabs performed testing. These are the conditions observed to be of most relevance to our customers. The dotted-line extensions to the solid-line segments represent extrapolated data and represent a wider range of conditions that may be encountered while operating the devices.

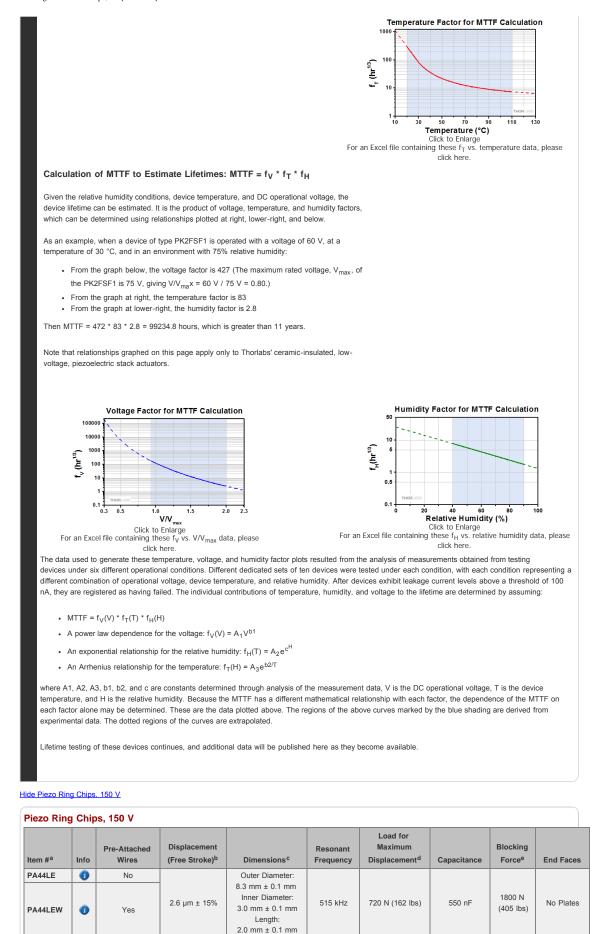

 All specifications are quoted at 25 °C, unless otherwise stated. We recommend using the MDT69xB or KP2101 Open-Loop Controllers with these piezos. Our MP2601 (≤75 V) or BPC30x Closed-Loop Controllers are also compatible. However, the actuators sold on this page do not contain a strain gauge and therefore will not provide positional feedback. If closed-loop feedback is needed. Delease consider our

Webpage Features Clicking this icon opens a window

containing item-specific specifications

PK4FYC2 or PZS001 actuators.

- · The "free stroke" displacement corresponds to no load.
- Dimensions are quoted for the chip. In the case of our wired products, the dimensions exclude the wire connection area.
- Displacement varies with loading. When used with this load, these chips achieve the maximum displacement, which is larger than the free stroke displacement.
- At Max Voltage

Part Number	Description	Price	Availability
PA44LE	Piezo Ring Chip, 150 V, 2.6 μm Displacement, 8.3 mm OD, 3.0 mm ID, 2.0 mm Long, Bare Electrodes	\$75.00 Volume Pricing Available	Today
PA44LEW	Piezo Ring Chip, 150 V, 2.6 $\mu m$ Displacement, 8.3 mm OD, 3.0 mm ID, 2.0 mm Long, Pre-Attached Wires	\$77.00 Volume Pricing Available	Today

# Hide Discrete Piezo Ring Stacks, 150 V

#### (\_\_. ..... zo Ring Stacks 150 V

Discrete P	Discrete Piezo Ring Stacks, 150 V								
Item # <sup>a</sup>	Info	Pre-Attached Wires	Displacement (Free Stroke) <sup>b</sup>	Dimensions <sup>c</sup>	Resonant Frequency	Load for Maximum Displacement <sup>d</sup>	Capacitance	Blocking Force <sup>e</sup>	End Face
PK44LA2P2	0	Yes	9.0 µm ± 15%	Outer Diameter: 8.3 mm ± 0.1 mm Inner Diameter: 3.0 mm ± 0.1 mm Length: 9.0 mm ± 0.1 mm	127 kHz	720 N (162 lbs)	2.0 µF	1800 N (405 lbs)	Flat Plate

· All specifications are quoted at 25 °C, unless otherwise stated. We recommend using the MDT69xB or KPZ101 Open-Loop Controllers with these piezos. Our MPZ601 (≤75 V) or BPC30x Closed-Loop Controllers are also compatible. However, the actuators sold on this page do not contain a strain gauge and therefore will not provide positional feedback. If closed-loop feedback is needed, please consider our PK4FYC2 or PZS001 actuators.

Webpage Features Clicking this icon opens a window containing item-specific specifications and mechanical drawings.

0

and mechanical drawings.

- · The "free stroke" displacement corresponds to no load.
- Dimensions are quoted for the chip. In the case of our wired products, the dimensions exclude the wire connection area, Kapton® tape, and wires.
- Displacement varies with loading. When used with this load, these chips achieve the maximum displacement, which is larger than the free stroke displacement. At Max Voltage

Part Number	Description	Price	Availability
PK44LA2P2	Discrete Piezo Ring Stack, 150 V, 8.3 µm Displacement, 9.7 mm OD, 3.0 mm ID, 9.0 mm Long, Two Flat End Plates	\$125.00 Volume Pricing Available	Today